
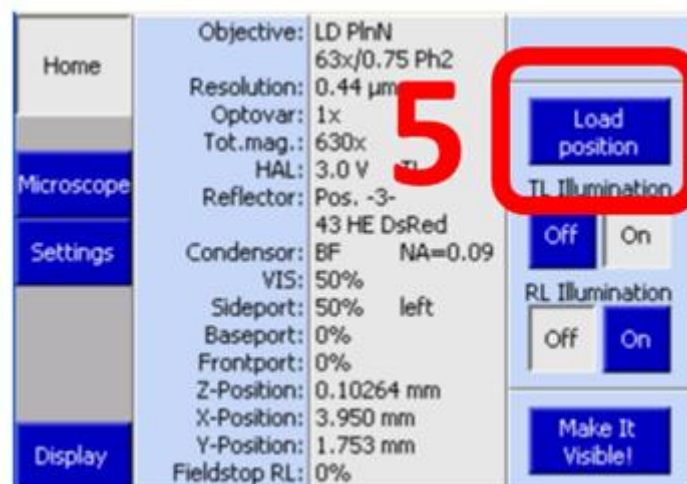


Startup

1. Turn the **Main Switch (1)** in ON position
2. Turn the **Laser key (2)** clockwise in vertical position (labeled "Laser")
3. Push the **Components** switch (3) in ON position
4. Power up the **PC (4)**. Login into the "user" account and start ZEN
5. Press the **Load position** button on the touch panel of the microscope

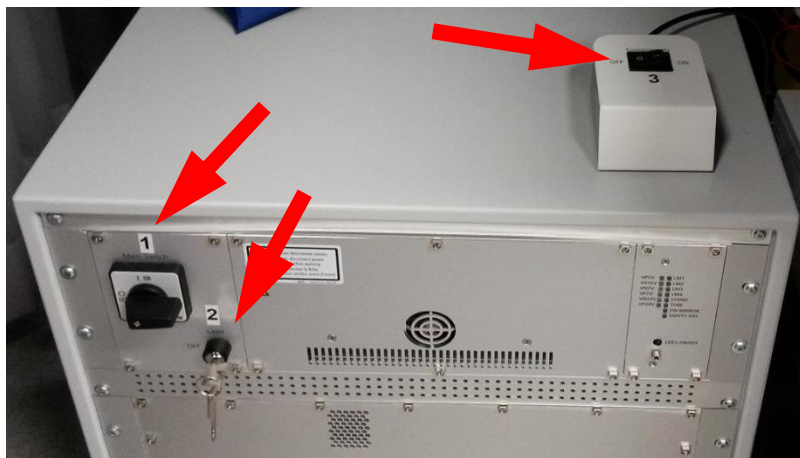

Shutdown

1. Close ZEN
2. Press the **Load position** button on the touch panel
3. Clean immersion objectives with lens cleaning paper

a. Leaving the system for the next user

4. Logout from the Windows session

b. Full shutdown (last user)

4. Shut down the **PC (4)**
5. Push the **Components** switch **(3)** in OFF position
6. Turn the **Laser key (2)** counter-clockwise in horizontal position (labeled "OFF")
7. Turn the **Main Switch (1)** in OFF position
8. Cover the microscope with the dust cover